

## **Application consideration**


### **RCB1514 : HOCE**

Please keep the following points in mind :

**Please DO NOT change medium right after thawing or next day of subculture, because these cells take time to attach to dishes.**

**Follicular formed aggregates will appear during cell culture(see the photo below). These cells can not be made into single cells by the treatment of trypsin-EDTA. Moreover, prolonged trypsin-EDTA treatment will cause some problems in growth after subculture. Therefore, when cells start to detach in sheet-like formations, please pipette up and down GENTLY and then plate cells to dishes.**

**Please subculture cells with high cell number density when cells grow in a polylayer manner even before cells reach to confluence.**

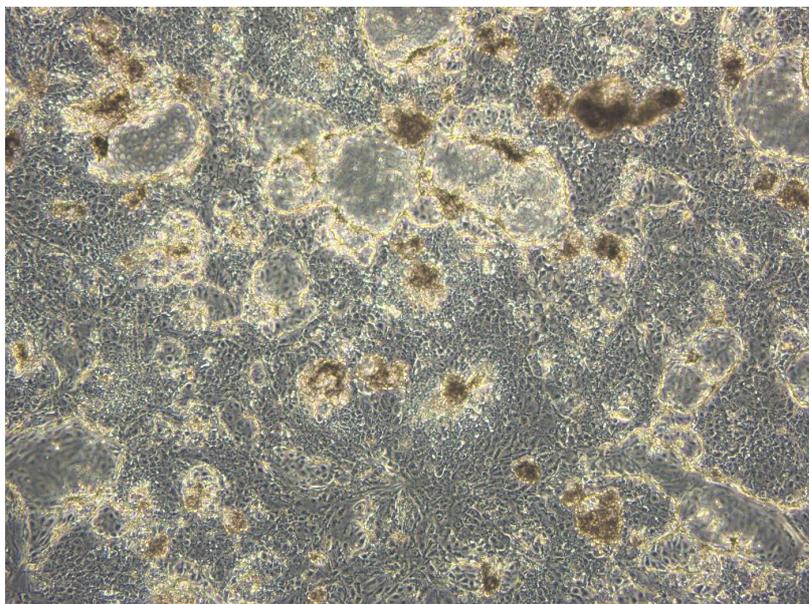

【Technical support】 e-mail : [cellqa.brc@riken.jp](mailto:cellqa.brc@riken.jp)